


Hydrolysis Unit B-411

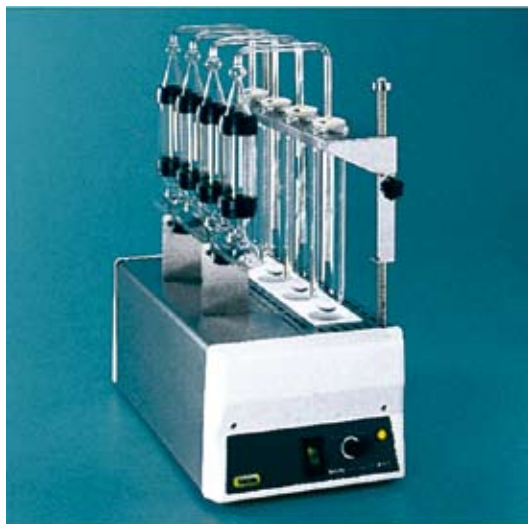

Hydrolysis



The ideal complement to the Extraction System

The Hydrolysis Unit B-411, offers the ideal complement to the Extraction System B-811. Collection of the entire fat with direct extraction is not always successful when doing fat determination according to Soxhlet. In such cases, hydrolysis before the extraction is necessary (Weibull-Stoldt). The Hydrolysis Unit B-411 provides problem-free acid digestion with minimal filtration time. Four determinations can be done simultaneously. The glass sample tubes are inserted directly in the Extraction System after drying.

Many different areas of use

The Hydrolysis Unit B-411, makes an important contribution to efficient fat determination, for example in food and animal feed or cosmetics. It is used in food and monitoring laboratories as well as in process monitoring.



Milk products



Meat



Cereal




Foodstuffs




Cosmetics

Hydrolysis

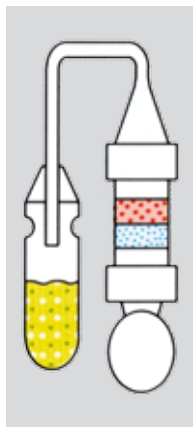

Filtration and rinsing

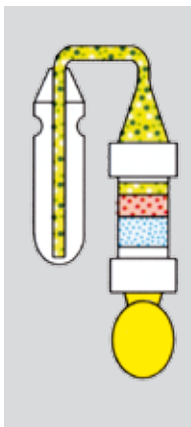

Clear-cut advantages

- Simultaneous hydrolysis of 4 samples
- Quick and efficient filtration
- User-friendly operation
- No odor nuisance
- AOAC and LMBG § 35 (G)

Extraction



The glass sample tube can be placed directly in the Extraction System B-811, for extraction after drying.

Delivery:

	Order No.
B-411 Hydrolysis Unit 230 V/50/60 Hz	37455
B-411 Hydrolysis Unit 120 V/50/60 Hz	37456
1 set of 4 digestion vessels 300 ml	37377
1 set of 4 glass sample tubes	37281
4 sample aspiration tubes	37380
1 set of 4 caps for digestion vessels	37463
1 set of 4 stoppers Ø 45	37725
1 inbus wrench 3.0 mm	00610

Accessory:

	Order No.
1 holder for 4 digestion vessels	02013
1 crucible stand (microwavable)	37462
1 water jet pump, plastic	02913

Technical data:

Number of heating points	4
Glass sample tubes (volume)	60 ml
Current consumption (220 V)	5 A
Power consumption	1100 W
Connected load	100–120 V, 220–240 V
Dimensions in mm (w x h x d)	275 x 600 x 570
Weight (net)	13.5 kg

BÜCHI Labortechnik AG
Postfach
9230 Flawil 1
Schweiz
T +41 71 394 63 63
F +41 71 394 65 65
buchi@buchi.com
www.buchi.com

BÜCHI Labortechnik GmbH
Postfach 10 03 51
45003 Essen
Deutschland
Freecall 0800 414 0414
T +49 201 747 490
F +49 201 237 082
deutschland@buchi.com
www.buechigmbh.de

BÜCHI Labortechnik GmbH
Branch Office Netherlands
Postbus 142
3340 AC Hendrik-Ido-Ambacht
The Netherlands
T +31 78 684 94 29
F +31 78 684 94 30
netherlands@buchi.com
www.buchi.nl

BÜCHI Italia s.r.l.
Centro Direzionale, Milano Fiori
Pal. A-4, Strada 4
20090 Assago (MI)
Italia
T +39 02 824 50 11
F +39 02 57 51 28 55
italia@buchi.com
www.buchi.it

BUCHI (Thailand) Ltd.,
77/175, Sin Sathon Tower,
39th FL., Unit F
Krungthongburi Rd.
Klongtonnai, Klongsan
Bangkok 10600
Thailand
T +66 2 862 08 51
F +66 2 862 08 54
bacc@buchi.com
www.buchi.com

BUCHI SMP
Services Private Ltd.
201, Magnum Opus
Shantinagar Industrial Area
Vakola, Santacruz (East)
Mumbai 400 055,
India
T +91 22 66 98 94 50 / 51
F +91 22 66 98 94 52
smplisp@vsnl.com
www.buchi.com

BUCHI Corporation
19 Lukens Drive, Suite 400
New Castle
Delaware 19720
USA
T +1 302 652 3000
F +1 302 652 8777
Toll Free: +1 877 692 8244
us-sales@buchi.com
www.mybuchi.com

BUCHI Hong Kong Ltd.
1810 Fortress Tower
250 King's Road
North Point, Hong Kong
China
T +852 2389 2772
F +852 2389 2774
china@buchi.com
www.buchi.com

BUCHI Shanghai Trading LLC
21/F Shanghai Industrial
Investment Building
18 Caoxi Bei Road
200030 Shanghai
China
T +86 21 6468 1888
F +86 21 6428 3890
china@buchi.com
www.buchi.com

BUCHI UK Ltd
5 Whitegate Business Centre
Jardine Way
(off) Broadway
Chadderton
Oldham OL9 9QL
United Kingdom
T +44 161 633 1000
F +44 161 633 1007
uk@buchi.com
www.buchi.co.uk

BUCHI Sarl
5, rue du Pont des Halles
Z.A. du Delta
94656 Rungis Cedex
France
T +33 1 56 70 62 50
F +33 1 46 86 00 31
france@buchi.com
www.buchi.fr

Nihon BUCHI K.K.
3F IMON Bldg.,
2-7-17 Ikenohata, Taito-ku,
Tokyo 110-0008
Japan
T +81 3 3821 4777
F +81 3 3821 4555
nihon@buchi.com
www.nihon-buchi.co.jp

We are represented by more than 100 distribution partners worldwide. Find your local representative at www.buchi.com

Quality in your hands